Dino-Lite

Digital Microscope

The Industry Standard

Dino-Lite Mobile User Manual

EN **Manual**
FR **Manuel d'utilisation**
DE **Bedienungsanleitung**
ES **Manual de usuario**
NL **Handleiding**

**Digital Video Recorder with
Dino-Lite digital microscope**

**AMK4012T-D15
AMK4012TL-D15
AMK4012ZT-D15**

www.dino-lite.eu

Redefine the microscope

AMK4012T/ZT/TL-D15

OPERATION MANUAL

EN	Manual	2
DE	Bedienungsanleitung	11
FR	Manuel d'utilisation	22
ES	Manual de usuario	34
NL	Handleiding	45

AMK4012T/ZT/TL-D15

OPERATION MANUAL

Digital Video Recorder with Dino-Lite digital microscope

Operating instructions

Before operating the unit, please read this manual thoroughly, and retain it for future reference.

Thank you for purchasing a Dino-Lite with mobile recorder.


Important safety information

- Avoid touching the lens with your finger to protect the product from electrostatic damage.
- Do not drop
- Keep dry
- Do not use it in high temperature, high humidity environment, 0~40°C is the best working temperature for this product
- Please insert or pull out the SD card or the Dino-Lite when the power is off. The product or the SD card could be damaged.

Please check www.dino-lite.eu for updated information or if you have any questions.

Basic Information

- **Display:** 5-inch HD screen (800X480pix).
- **Channel Switch:** Supports manual & automatic switch between 8 channels.
- **Recording:** Supports seamless cycle and continuous video from the Dino-Lite digital microscope
- **Photo:** Supports function of taking pictures either from the Dino-Lite digital microscope
- **AV IN:** TV-IN [NTSC/PAL].
- **AV OUT:** NTSC/PAL].
- **Battery :** Built-in 3200mAh high capacity Li-battery, continuous record over 10 hours.
- **Speaker:** Built-in speaker of 1W.
- **Storage:** Supports up to 32GB TF card, 4GB SD card supplied.
- **Language:** Supports English/ French/German/Spanish/Dutch.
- **Power-saving Mode:**Automatic shutdown, automatic screen off.
- **Firmware update:** Supports firmware update from SD card.

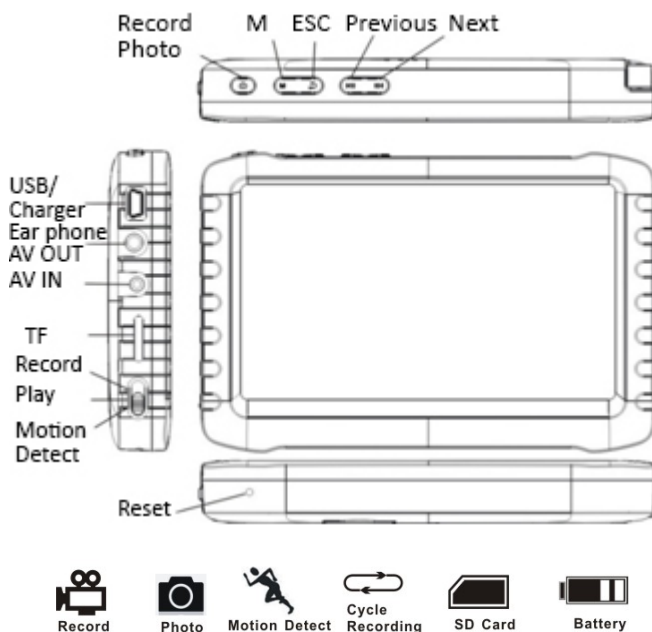Connecting the Dino-Lite digital microscope

With the recorder power set to off, connect the plug on the Dino-Lite to the AV-IN connector on the recorder.

Note: The AMK4012 Dino-Lite is not supported for use on other devices with 2,5 mm AV connector.

How to Charge

Please make sure to charge the battery fully before first usage. Only use the charger supplied with the product or charge on a PC USB port.

Charge 4~6 hours for a full charge. Do not unplug the charger until the battery status is full. For the first time charge, please charge for more than 6 but less than 10 hours. A new battery will reach best performance after a few times full charge and discharge. If the recorder is not used for a longer period, please recharge every 2-3 months.

Battery capacity symbols:

Basic Operation

1. Power On/Off: Press “**M**” to power on, DVR is in working state; press “**M**” for several seconds to power it off.

2. Working mode switch (switch located at the left corner):

When the switch is on “**📷**”, DVR is in state of receiving/record/photo ;

When the switch is on “**▶**”, DVR is in playback;

When the switch is on “**🚗**”, DVR is in the state of motion detection.

3. External Storage card

This DVR can only record photo or video if a storage card has been plugged in. The screen will show symbol of SD card & its free capacity.

4. Working as a disk

Connect DVR to computer and power it on, it will work as a disk. Open the removable disk in the computer, you'll find the videos and photos in “RECORD” and “PHOTO” files under “DCIM” directory.

Record & Photo

Power DVR on, and select working mode “” for recording/photo mode.

1. Photo: In “” mode, press “” once, symbol “” flashes once, and one picture is made.

2. Record: In “” mode, press “” for 3 seconds, the DVR will start to record and the screen shows “REC”. Press “” again to pause, screen shows “||”; press “” one more time, to resume recording. Press “” to end, it will save automatically and exit. If you do not press “”, the system will automatically save video in segments until the storage card is full. If “Overwrite” is turned on and the card is full, it will overwrite the earliest recording and save the latest one.

3. Mute: In “” mode, press “” to enter state of mute. Screen displays “MUTE”. Speaker gives no sound; press “” again to recover sound.

Note: Mute setting will not affect earphone volume.

Motion Detect

1. Power DVR on, and select working mode to “”, for motion detect mode. Press “” to exit and go back to system/function menu.


2. In “” mode, press “” for 3 seconds to enter into motion detect recording. When moving objects appear, system will automatically start to record, and screen will show “REC”. After 10 seconds without moving objects, system will automatically end recording and save to exit, waiting for next motion. To end motion detect recording, press “”, it will save automatically and exit.

3. If you do not press “”, the system will automatically save video in segments until the storage card is full. If “Overwrite” is turned on and the card is full, it will overwrite the earliest recording and save the latest one.

Play Video/Photo

When the working mode is set to "⏮" you can playback video & photo. Press "⏮"/"⏭" to select "Photo" or "Video". Press "⏮" to exit and go back to main menu.

1. Play Video: When you select "Video", press "📷" to enter into date folder. Press "⏮"/"⏭" to select the date, and press "📷" to enter into the list of video files. Press "⏮"/"⏭" to select, and press "📷" to play video. Press "⏮"/"⏭" to adjust volume. Press "⏮" long time to fast forward; press "⏭" long time to rewind; press "📷" to pause; press "📷" again to release pause; press "⏮" to exit.

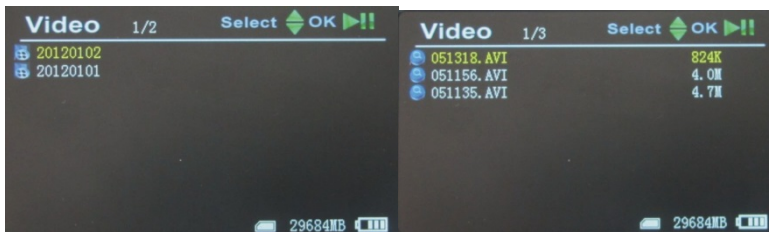2. Play Photos: Select "Photo" and press "📷" to enter into photo folder. Press "⏮"/"⏭" to select pictures, then press "📷" to view. Press "⏮" to view next one; press "⏭" to view previous one. Press "⏮" to exit.

3. Delete Photo & Video: In list mode, press "⏮"/"⏭" to select files. Press "M" to get the delete dialogue. Press "⏮"/"⏭" to select "All"/"Current"/"Cancel". Press "📷" to confirm.

Function Settings

Power DVR on, working mode switched to "📷" / "⏮", press "⏮" to enter into function and system setting menu. Press "⏮"/"⏭" to select "Function", then press "📷" to enter into function menu.

1. Recorder Resolution: Press "⏮"/"⏭" to select "Recorder Resolution" and press "📷" to enter. Press "⏮"/"⏭" to select "720X576" / "720X480". Press "📷" to confirm and exit.

2. Frame Rate: Press “”/“” to select “Frame Rate” and press “” to enter. Press “”/“” to select “30 FPS”/“25 FPS”, and press “” to confirm & exit.

3. Video Segment Sizes: Press “”/“” to select “Video Segment Sizes”, then press “” to enter. Press “”/“” to select “10 MIN”/“30MIN”/“60 MIN”. Press “” to confirm & exit.

4. Overwrite: Press “”/“” to select “Overwrite”, and press “” to enter. Press “”/“” to select “Off”/ “On”, and press “” to confirm & exit. If “Overwrite” is on, DVR will cycle record, and save the latest video.

5. Time Stamp: Press “”/“” to select “Time Stamp”, and press “” to enter. Press “”/“” to select “Off”/ “On”, and press “” to confirm & exit. If it’s on, video will be saved with synchronous time.

6. Information Display: Press “”/“” to select “Information Display”, and press “” to enter. Press “”/“” to select “Off”/ “On”, and press “” to confirm & exit. If it’s on, screen will display information of working frequency, channel, audio state, date & time, etc. If you select “Off”, no information will be displayed.

7. Video Setting: Press “”/“” to select “Video Setting”, and press “” to enter. Press “”/“” to select “Brightness” or “Contrast”; then press “” to enter; then press “”/“” to adjust brightness or contrast. Press “” to exit

System Settings

Power DVR on, working mode switched to “” / “”, press “” to enter into function & system setting. Press “”/“” to select “System”; then press “” to enter into system menu.

1. Date time: Press “”/“” to select “Date Time”, and press “” to enter. Press “” again to move the cursor, and press “”/“” to adjust date & time. Press “” to exit.

2. Language: Press “”/“” to select “Language”, and press “” to enter. Press “”/“” to choose your language. And press “” to confirm and exit.

3. AV Out: Press “”/“” to choose “AV Out”, and press “” to enter. Press “”/“” to choose “PAL/NTSC”, press “” to confirm video displayed on TV screen. Pull out AV cable, video will show on DVR screen.

4. Format: To format the SD card, press “”/“” to select “Format”, and press “” to enter. Press “”/“” to choose “Yes”. Press “” to confirm & exit.

5. Power Save: Press “”/“” to select “Power Save”, and press “” to enter:

a) Auto LCD Off: Press “”/“” to select “Auto LCD Off”, and press “” to enter. Press “”/“” to select “1 MIN”/“5 MIN”/“Off”, and press “” to confirm & exit. If “Auto LCD Off” is on, the LCD screen will automatically shut off when the set time is due. Press any key to recover screen display.

b) Auto Power Off: Press “”/“” to select “Auto Power Off” and press “” to enter. Press “”/“” to select “5 MIN”/“10 MIN”/“Off”, and press “” to confirm & exit. If you turn on “Auto Power Off”, the DVR will automatically shut off when the set time is due.

6. Firmware Update: Press “”/“” to select “Firmware Update”, and press “” to enter. Press “”/“” to select “Yes”, and press “” to start update. Please remember to save the update file “zburnT_C11.run.elf” at the root directory of the SD card.

7. Default Setup: Press “”/“” to select “Default Setup” and press “” to enter. Press “”/“” to select “Yes” and press “” to confirm. Press “” to exit.

8. Memory Info: Press “”/“” to select “Memory Info”, and press “” to view capacity of SD card. Press “” to exit.

9. Version: Press “”/“” to select “Version” and press “” to view version information. Press “” to exit.

10. Display Mode: Press “”/“” to choose “Display Mode”. Press “”/“” to choose “16:9/4:3”. Then press “” to confirm. Press “” to exit.

Troubleshooting

1. The recorder does not power on: maybe the battery is empty, try to recharge.
2. The Dino-Lite is connected but there is no image: check if the plug is in the AV-in port and properly inserted. If this does not work, try to change the settings or return to Default Setup.
3. System halted: Press the RESET key to turn on player again.
4. File cannot play: Maybe the format of video file is not supported by this DVR, please convert it into the supported format.
5. The text is in messy code: make sure the correct language is selected.
6. It can't upload / download files or can't find the disk in PC: Replace USB cable/USB port.
7. It cannot save a file: there is no SD card in theDVR, please insert SD card; or SD card memory is full, please delete some files to release memory; or there is damaged area in memory, please format it.

Package Content

The following items are in the package.

Dino-Lite AMK digital microscope	1x
MSD15 recorder	1x
Clip	1x
User manual	1x
USB Cable	1x
AV Cable	1x
AC adapter(with EU/UK/US plug)	1x
4GB Micro SD Card	1x

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Warranty

The Dino-Lite product is delivered with a two year warranty from the date of purchase by the end customer. The MS-D15 recorder is delivered with a one year warranty from the date of purchase by the end customer. Please note that accessories (i.e. detachable caps/extensions) are not covered under warranty. For warranty issues please contact the reseller or store where you bought the product.

The content of this manual has been prepared with care, but Dino-Lite Europe does not warrant for any mistakes or omission. Dino-Lite Europe is not responsible for loss of personal data, productivity or otherwise resulting from wrong operation or product failure. The software, hardware and this manual may be modified without further notice.

Support

If you have any problem or issue with your Dino-Lite or the MS-D15 recorder, please contact your reseller or check the Dino-Lite Europe support pages at www.dino-lite.eu.

Dino-Lite Europe/IDCP BV
The Netherlands
Tel. +31 20 6186322
Fax. +31 20 6189692
e-mail: info@dino-lite.eu
web: www.dino-lite.eu

AnMo Electronics Corporation and its authorized representative in the EU, IDCP BV, declare that the apparatus Dino-Lite comply with the essential requirements and other relevant provisions of Directive 1999/5/EC. The original Declaration of Conformity can be found at www.dino-lite.eu/doc/dino-lite.pdf.

IDCP BV declares that the MS-D15 recorder complies with the essential requirements and other relevant provisions of Directive 1999/5/EC. The original Declaration of Conformity can be found at www.dino-lite.eu/doc/msd15.pdf.

2014 © AnMo/IDCP BV/Dino-Lite Europe.

No part of this publication may be reproduced or distributed unless for personal use in conjunction with usage of Dino-Lite digital microscopes.