

# Dino-Lite Pro AMK412-C200

Dino-Lite Pro AMK412-C200 at a glance



AMK-412

## Accessory



C200



RCA Cable










Battery

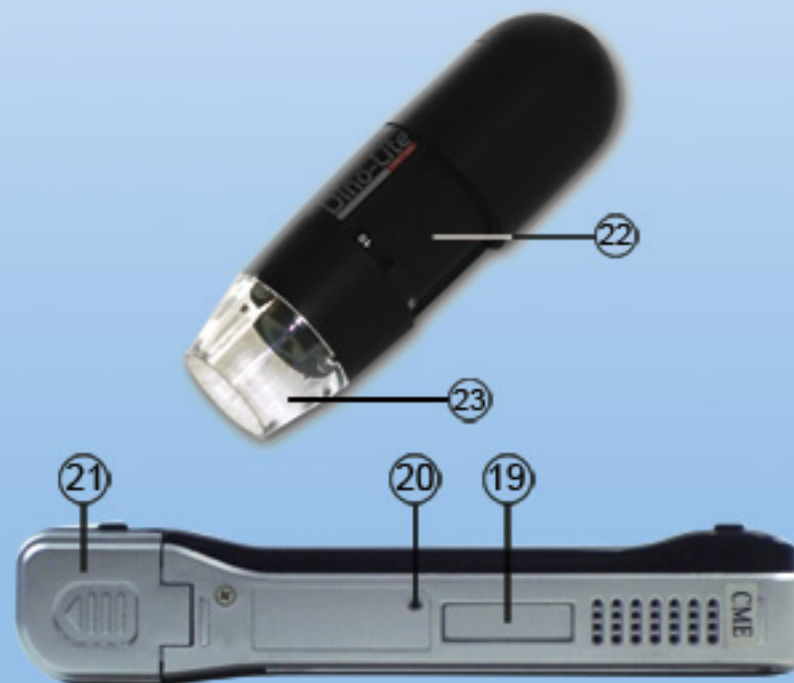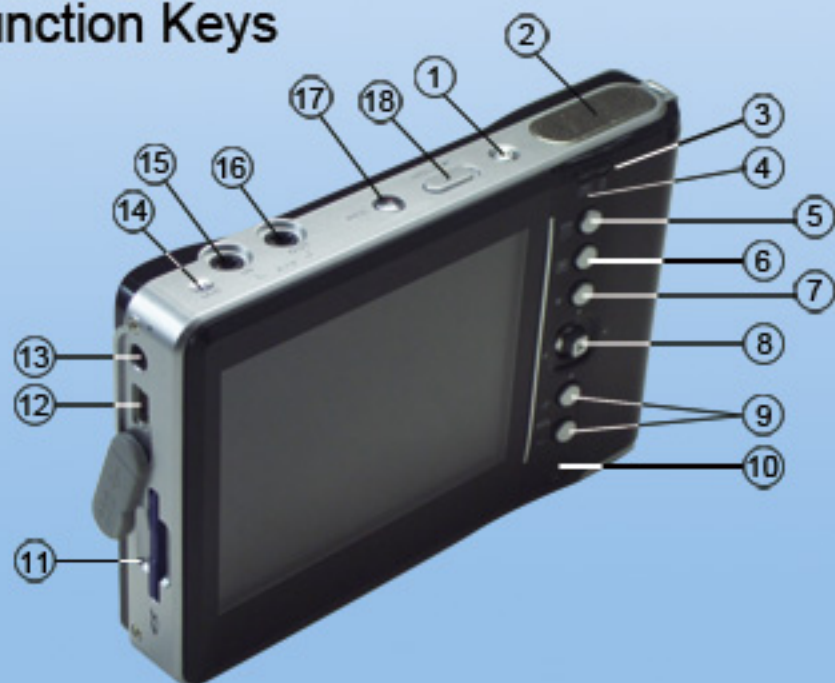

Battery Charger

## Function Keys



1. Power Switch	9. Volume Control 	17. A/V Recording
2. Speakers	10. Remote Control Receiving Interface	18. Key Lock
3. Charging Status Indicator	11. SD Card Slot	19. Expanding Slot Connector
4. Recording Status Indicator	12. USB Slot	20. Reset Key
5. Setup 	13. Battery Charging Slot	21. Battery Cover
6. Snap 	14. Microphone	22. Focus & Magnification Adjustment Knob
7. Stop 	15. AV-Input	23. LED Lighting & Cover
8. 5-way Navigation Keys 	16. AV-Output / Earphone jack	

# Dino-Lite Pro AMK412-C200

## Charging the Battery of C200

For the first time user, please charge the battery of C200 for more than 8 hours by connecting C200 to the adapter and then connect the adapter to the power outlet. The charging can also be performed by connecting the C200 to a computer with an USB cable.

### Notice:

The C200 may not function when the battery is empty, even the adapter is connected.

To extend the battery's life, please unplug the adapter immediately when the charging status indicator LED is vanished.


## Starting up the AMK412-C200


1. Connect Dino-Lite AMK412 to AV-Input of C200 before power on the C200.
2. Make sure the SD card is inserted in the SD card slot of C200.
3. Turn on the C200 by pressing the power switch. The LED of AMK412 should be then turned on, and the following screen will be displayed initially for 2 seconds and the screen will then display the realtime video transmitted from the AMK412.


## Taking Picture




Taking picture can be done by simply pressing . The taken picture will be stored in the SD card.

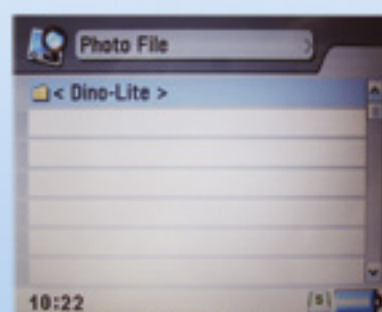

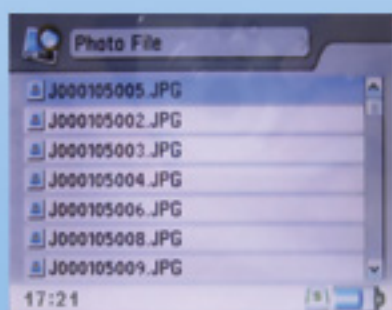

To review the picture, pressing  for one time to enter the Photo File. A folder of <Dino-Lite> can then be found.




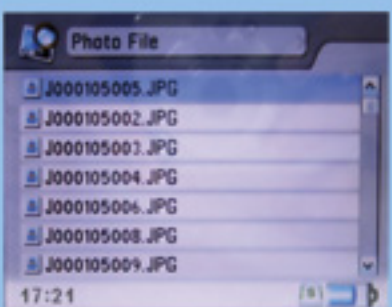
Go into the <Dino-Lite> folder by pressing , and a list of picture files will be shown as below.




Scrolling the Photo File List by using  to the file to be selected, then press  and the picture will be displayed.




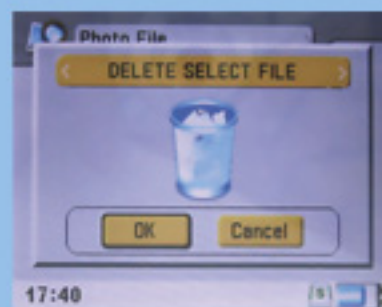
The picture can be 90-degree rotated by pressing .



Return to the Photo File List or the preview screen by pressing .





Delete the picture in the Photo File List can be done by pressing  and the following screen will be displayed.




Delete single picture or all pictures by selecting One or All in the delete option.

## Taking Video




Taking video can be done by simply pressing AV Recording button, and press again to pause recording. Stop or return from the recording by pressing . Voice is able to be recorded during the video is taking with C200' s built-in microphone. The taken video will be stored in the SD card.

To review the video, pressing  for **two times** to enter the Video File. A folder of <Dino-Lite> can then be found.




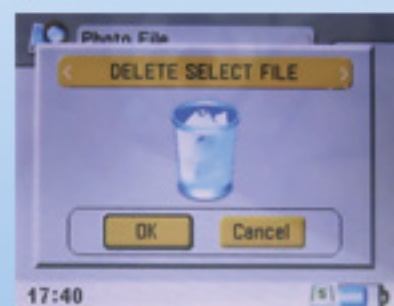
Go into the <Dino-Lite> folder by pressing , and a list of video files will be shown as below.



Scrolling the Video File List by using [: up/down] to the file to be selected, then pressing  and the video will be displayed. Return to the Video File List or the preview screen by pressing .




Delete the video in the Video File List can be done by pressing  and the following screen will be displayed.





Delete single video or all video by selecting One or All in the delete option.

## Setup

The setup provides functions including setting of time, TV format, display, language, and also format of SD card, backup from SD to harddisk if a harddisk in equipped, and firmware upgrade etc.

By pressing  for three times in the preview mode, a setup screen will then be displayed as following.



Use  to select the setup function and change it' s setting. Return from Setup by pressing  or untouch any key for more than 3 seconds.